



P.O. Box 2307
Westerville, OH 43086
(614) 895-2837
www.Mindwaretech.com

**!! DO NOT INSERT THE USB KEY INTO THE USB PORT UNTIL
INSTRUCTED TO DO SO !!**

Installation Instructions

To install the MindWare applications it is important that you follow these instructions in the correct sequence. There are three steps to the procedure that include installing Nidaq, installing the application, and installing the USB protection key. This is a quick process that should have you up and running in 15 minutes.

Step 1: Installing Nidaq from the provided cd

If Nidaq is already installed on your pc, you do not need to reinstall it. This step is only for those users who are installing it for the first time.

Insert the Nidaq cd into your cd-rom drive. It is set to autorun and should launch automatically. If for some reason it doesn't automatically launch, you can simply run the setup.exe program on the cd. Follow the default responses as the program executes. A reboot will be required to complete the installation.

Step 2: Installing the MindWare application

Insert the MindWare application cd into your cd-rom drive. As with the Nidaq installation, it is set to autorun and should launch automatically. It is recommended that you keep the default parameters as the application installs.

The MindWare applications require that the provided USB key be present to run. ***Important: DO NOT INSERT THE USB KEY INTO THE USB PORT UNTIL AFTER THIS STEP OF THE INSTALLATION FINISHES.*** The required hardware drivers for the key will be installed so that the New Hardware Found Wizard can load them when required.

Step 3: Inserting the USB Key (Not required for Demo Version)

Now is the appropriate time to insert the provided USB key into one of the USB ports on the pc. Once the operating system detects the device it will launch the New Hardware Found Wizard to finish the installation. Once this has completed, you are ready to run your MindWare application.