

## RELEASE NOTES: LXSoftware v 11.10.3

Released: October 29, 2024

### Minor Updates

- Remove PolyScore
- Modify RLE constraints to natural log of 1.2 – 1.6 in both positive and negative directions
- Replace the obsolete Functionality Check hardware manual with document describing how to perform a Field Functionality Check

### Bug Fixes

- Resolve issue that when scoresheet was configured to show only one pneumo sensor the pneumo score was not added to the subtotal

## RELEASE NOTES: LXSoftware v 11.10.2

Released: August 21, 2024

### Bug Fixes

- Resolve issue with NCCA import that Automatic EDA data would be incorrectly imported as Manual EDA data

## RELEASE NOTES: LXSoftware v 11.10.1

Released: August 15, 2024

### Minor Updates

- If not already present, automatically install Microsoft Visual C++ 2015-2022 Redistributable Package
- For NCCA export, stop adding “R” or “C”, if missing, to Relevant or Comparison question labels, respectively
- Added the Import and Export buttons back to the Question Manager

## RELEASE NOTES: LXSoftware v 11.10.x

### Bug Fixes

- Resolve issue that “Documents and Attachments” menu item did not bring up the expected window
- Resolve issue that duplicate PFs (ending in ~1, ~2, etc.) were being created when opening PF using file explorer or folder browser

## RELEASE NOTES: LXSoftware v 11.10.0

Released: August 6, 2024

### Major Updates

- Update Build tools to MVC142
- Update to C++17
- Downgrade Video Capture library to version used in 11.8.6.399; this resolves significant issues that were seen during video and audio recording
- Prevent installing over existing LXSoftware install
  - If user attempts to install over existing LXSoftware, displays message box with instructions

### Minor Updates

- ESS
  - Correctly compute the number of control questions on old PFs
  - Ensure that Prior Odds appear in the ESS report and in the PDF
- Masseter Headset
  - Emplace UI to allow selection of the headset under Preferences->Admin
- Allow for registry configuration to prevent usage of FTP or scripting
- Remove support for the Dari language
- Change default Activity Seat gain from 5.0 to 1.0
- Stop resizing Recorder application when selecting a different camera
- Add the OSS-3 manual to the Polygraph References library
- Remove “Transfer Files via FTP” from the Tools menu
- Removed the Import and Export buttons from the Question Manager

### Bug Fixes

- ESS
  - Resolve chart selection error
  - Fix issue that extra text would sometimes appear in the ESS Report title
- Masseter Headset

## **RELEASE NOTES: LXSoftware v 11.10.x**

- Resolve class of merge errors due to timing differences
- Fix setting that turns off warning that appears when creating new series
- Fix minor issues with text appearance when adjusting gain
- Resolve issue that could cause LXSoftware to freeze
- Fix the scoring artifact character in Spanish and Russian versions